

PLANER Shipslog Software Download Instructions

1 Introduction

The **ShipsLog**[™] is a temperature monitoring system for samples transported in liquid nitrogen vapour shippers. It provides traceability for the temperature of the samples and can provide alarms at predefined setpoints during the transit period.

The **ShipsLog** monitor is supplied either fitted to the top of the vapour shipper's cover or with a flying lead for positioning within the sample area.

Samples may be transported in the knowledge that on arrival at the destination, the recipient can quickly check the LED display and confirm whether the temperature experienced by the samples has deviated outside of the predefined limits. The full temperature history can also be downloaded to confirm the exact temperature pattern throughout the journey.

The logger is user-programmable via PC software. The logging process can be started either by using the PC software or by using the magnetic swipe key provided. By using the PC software, the data can be analysed by both sender and receiver.

2 Installing the software

In order to use the unit, the **ShipsLog for Windows 32** software must first be installed. This software can be downloaded from the Planer website at <http://www.planer.com>. From the website, follow the link to **Support > Downloads > Products > ShipsLog** section to locate the download link for the **Shipslog** software.

When you have downloaded the file, unzip the files to a suitable temporary folder on your PC. To install the software, execute the file **Setup.exe**, located in the folder where you unzipped the files and then follow the on-screen instructions.

When the installation has completed, launch the **ShipsLog for Windows 32** application and then select **Help | ShipsLog Operation** from the main menu.